



Priorclave details - model no PS/RSC/SH230.

SH230 RSC Priorclave

Built to a compact design allowing them to fit through a standard size door these autoclaves are available in 230, 285 & 350 litre capacities with other sizes available on request. Castor mounted for ease of positioning, they are available either electrically or direct steam heated and feature a simple, two hand wheel hinged door closure. All models are fitted with two non-tipping shelves and a range of stainless steel baskets, waste containers and loading trolleys are available to suit these autoclaves.



Rectangular Section Priorclaves.

For those occasions when only 'square' will do! Perfect for high volume throughput as well as large or bulky items. These autoclaves combine a huge internal capacity with a small floor area and are ideally suited for use in a wide range of sterilising applications such as media preparation, liquids and diluent, waste, glassware, instruments, apparatus and health care.

They are extensively used within many industries, including food, drink, dairy, pharmaceutical, agricultural, education, healthcare, research establishments and industry.

Features include:

- Capacities from 230 to 700 Litres and above. Intermediate sizes available to suit individual locations where required.
- Operation up to 140C - 2.5 Bar
- Large rectangular chamber within a compact frame
- Steam generated within the chamber (electrically heated models) - Direct steam heated models available
- Complete with thermocouple entry and test ports
- Forced air cooling as standard helping to reduce cycle times
- Automatic timed freesteaming as standard for improved air removal and sterilising performance
- Delayed start and media warming function as standard for ready to pour media at the start of the day
- A large selection of options and accessories is available to help to get the most from your Priorclave autoclave
- Supplied with two full width non-tipping stainless steel shelves.
- Stainless steel baskets, waste containers and loading trolleys available

All models are fitted with the **TACTROL** microprocessor control system. After setting your required operating parameters press the start button for a fully automatic cycle. The control panel shows cycle status with digital readouts of temperature and time that can be viewed simultaneously. Temperature and time settings are fully adjustable within the operating range of the autoclave.

Power door and double door (pass-through) versions are also available along with full steam/water jackets if required.

Safety

Pressure vessels are fully **CE** marked and insurance approved to the European PED 97/23/EC. The autoclaves fully conform to BS2646-1993, BS5500 Cat.3, and EMC directives, amongst others.

All autoclave doors are fully interlocked for safety with a special safety handwheel mechanism, which cracks the gasket seal allowing any residual pressure to safely escape. A thermal cooling lock, preventing the door from being opened until the contents have cooled to a safe temperature is also fitted along with a pressure interlock.

A safety blow down valve and a low water level sensor to guard against boil-dry conditions (electrically heated models) afford further protection.

Epoxy coated panels and frame members are treated with the unique **Biocote** anti-bacterial agent, which is effective against all bacteria and fungi including MRSA.

Installation Requirements

Electrical: 3 Phase – 400V, 21kW, 30A (230, 285 & 350L Models), 400V,32.5kW, 60A (450L Model), 400V, 42kW, 60A (700L Model)

Steam: (Direct steam heated models only): - Dry Steam reduced to 40psi. ½ in BSP.

Drainage: - The single 1 in vent outlet pipe can be directed to a nearby drain. For make safe processing the discharge should be directed to a sealed discharge system; the system should lead by direct connection to a building drain or catchment tank.

For more specific advice about your particular installation requirements please contact our Technical Department or download our data sheet outlining the installation requirements for Priorclaves or any other type of Laboratory autoclave at: <http://www.priorclave.co.uk/downloads/installation.PDF>

Rectangular Section Priorclave autoclaves have a 12 months parts and labour warranty (parts only outside the UK) and a 10 year or 8,000 cycle pressure vessel warranty.

Technical Specifications.

Max. Capacity: 232L

Internal Diameter: -mm

Load height: -mm

Dimensions (mm)	Width	Depth	Height	Packed Dimensions (mm)	Width	Depth	Height
Internal	580	600	675	UK Delivery	770	1000	1660
External	760	950	1635	Delivery by Sea	-	-	-
Installed	960	1635	1150	Delivery by Air	-	-	-

Weights (Kg)	
Gross	500
Packed (Sea)	600
Packed (Air)	-

NB: Weights and dimensions shown are for the Autoclave only, unloaded and without water (if applicable). These figures do not include any options and accessories listed below which may alter external dimensions and weights. For full details please contact Priorclave or your local agent.
